


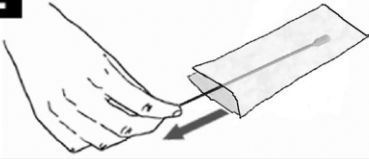

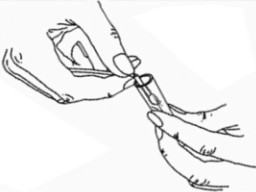

Instructions for Use: Isohelix BuccalFix RNA 0.5ml collection tubes for use with polyester breakable Buccal Swabs

REF RFSS-S1-05



Intended for the retrieval and stabilisation of buccal RNA. Single use only.
Store at room temperature. Use only if swab wrapper remains intact.

Precautions: Do not ingest. If the stabilising solution comes into contact with skin or eyes, wash with plenty of water, Cap may be a choking hazard to small children. MSDS available at www.isohelix.com.

<p>Take your RNA sample at least one hour after eating, drinking or cleaning your teeth. For best results, rinse mouth with water immediately prior to sampling.</p>	
1	 <p>Pull open the package from one end.</p>
2	 <p>Remove the swab from the wrapper, taking care not to touch the swab head with fingers.</p>
3	 <p>Insert the swab into the mouth and rub firmly against the inside of the cheeks or underneath the lips. Rub for at least 1 minute to maximise sample collection. Apply firm, reasonable pressure when swabbing.</p>
4	 <p>Remove cap from collection tube, place swab into tube and remove swab shaft either by snapping the swab head off, or by cutting off the swab head.</p>
5	 <p>Replace the tube cap, making sure the tube is securely sealed. Shake well to ensure that swab mixes with the collection buffer.</p>

RNA stability on swabs stored in BuccalFix RNA tubes at room temperature is maintained for two weeks.

FOR RESEARCH USE ONLY

Isohelix is a division of Cell Projects

For swab or DNA isolation queries email: info@isohelix.com www.isohelix.com
Molecular Biology Solutions www.cellprojects.com

Version August 2020