


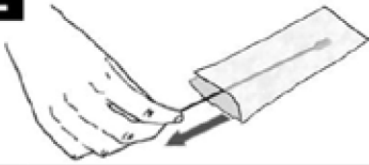



## Instructions for use of Isohelix RD-01 Buccal Swabs with RapiDri Pouch

REF RD-01

CE UK MD

Intended for the retrieval of buccal cells. Single use only.

Store at room temperature. Use only if swab wrapper remains intact.

<p><b>Take your DNA sample at least 1 hour after eating, drinking or cleaning your teeth.</b> <b>For best results, rinse mouth with water immediately prior to sampling.</b></p>	
<p><b>1</b></p> 	<p>Pull open the package from one end.</p>
<p><b>2</b></p> 	<p>Remove the swab from the tube, taking care not to touch the white swab head with your fingers</p>
<p><b>3</b></p> 	<p>Insert the swab into your mouth and rub firmly against the inside of your cheek or underneath lower and upper lip. For standard DNA collection rub for <b>1 minute</b> and in all cases rub for a minimum of 20 seconds. <b>Important – use reasonable, firm and solid pressure</b></p>
<p><b>4</b></p> 	<p>Place the swab head first into the RapiDri pouch so that the head sits within the drying area.</p>
<p><b>5</b></p> 	<p>Fold the RapiDri pouch at the fold line, with the clear side inwards as shown, and use the barcoded sticker to seal.</p>

It is the sole responsibility of the Laboratory or Professional User to verify and correctly validate the use, technically and legally, of these DNA Buccal swabs for their intended device or process.  
In case of any serious issues contact the manufacturer or local competent authorities.