


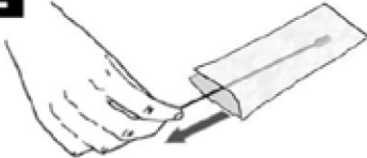



Instructions for use of Isohelix RD-01 Buccal Swabs with RapiDri Pouch

REF RD-01



Intended for the retrieval of buccal cells. Single use only.

Store at room temperature. Use only if swab wrapper remains intact.

<p>Take your DNA sample at least 1 hour after eating, drinking or cleaning your teeth. For best results, rinse mouth with water immediately prior to sampling.</p>	
<p>1</p> 	<p>Pull open the package from one end.</p>
<p>2</p> 	<p>Remove the swab from the tube, taking care not to touch the white swab head with your fingers</p>
<p>3</p> 	<p>Insert the swab into your mouth and rub firmly against the inside of your cheek or underneath lower and upper lip. For standard DNA collection rub for 1 minute and in all cases rub for a minimum of 20 seconds. Important – use reasonable, firm and solid pressure</p>
<p>4</p> 	<p>Place the swab head first into the RapiDri pouch so that the head sits within the drying area.</p>
<p>5</p> 	<p>Fold the RapiDri pouch at the fold line, with the clear side inwards as shown, and use the barcoded sticker to seal.</p>

It is the sole responsibility of the Laboratory or Professional User to verify and correctly validate the use, technically and legally, of these DNA Buccal swabs for their intended device or process.