


Do not eat, drink, smoke, brush your teeth or chew gum for 30 minutes before use.
Check fluid level is at 1ml line before opening. Inspect swabs for any damage before use.

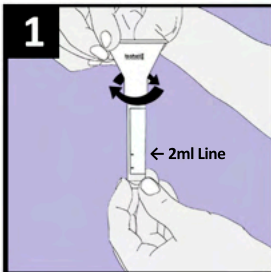

1. Remove the cap from the tube.
Keep cap for later use.
Screw collection funnel on to the tube.
Remove foam swabs from wrappers.

Do not leave donor unattended during sample collection.




2. Ensure donor is sitting in an upright position.
Gently insert swab into mouth and move around the front area of the mouth and under the tongue until swab is completely wet.

DO NOT - Chew, pull swab head, or use swab towards the back of the mouth.



3. Remove swab from mouth, and squeeze swab head against the side of the funnel using a rotating motion to dispense saliva from the sponge.

Once saliva is dispensed, repeat steps 2 and 3 multiple times until fluid in tube **reaches the 2ml line**.
If swab shows signs of wear and tear, use the second swab.

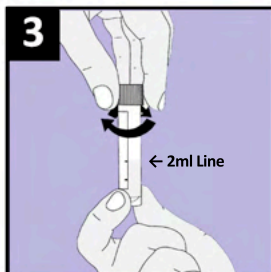

4. Unscrew the collection funnel and replace the tube cap tightly. Discard the funnel.

Shake the collection tube several times to mix the saliva with the solution.

STORE AT ROOM TEMPERATURE AFTER COLLECTION

Precautions: Do not swallow. If stabilising solution comes into contact with skin or eyes, wash with plenty of water. Tube cap and swabs may be a choking hazard to small children. MSDS available at www.isohelex.com.

In case of any serious issues contact the manufacturer or local competent authorities.